

CARSON[®] MicroBrite™ Plus

MM-350

60-120x LED Lighted Pocket Microscope with Adapter Clip

Congratulations on selecting your new MicroBrite™ Plus Microscope with Smartphone Adapter Clip!
In order to achieve optimum performance, please follow instructions for proper use and care.



PRODUCT SPECIFICATION


Magnification:	60x - 120x	Weight:	0.13 lbs (excluding battery)	Battery:	1 AA (not included)
Exit Pupil Diameter:	10mm - 14mm	Dimensions:	2.3" x 1.2" x 3.7"	Accessories:	MM-310 Clip
Optical Zoom:	2x	Illumination:	LED		

INSTALLING BATTERY

1. Slide open battery door.



2. Insert 1 AA battery with the correct polarity and replace battery door.

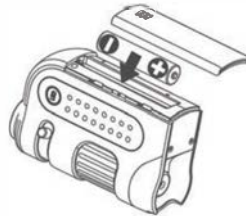


WARNING: Use only alkaline battery. Keep away from fire or excessive heat. Inspect battery for corrosion or leakage. Insert in direction indicated. Rechargeable battery not recommended.

USING MICROSCOPE

1. Place viewing object directly below Objective Lens.



2. Press LED Button to turn ON/OFF.

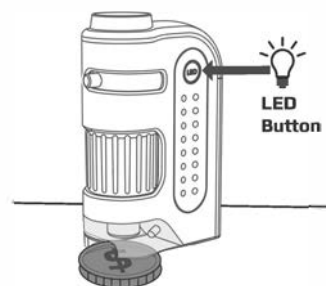

3. Set Zooming Dial to lowest magnification and rotate Focusing Ring slowly until image is clear. Adjust zoom and refocus as needed.



Note: To work properly, microscope must be placed flush to viewing object.

USING ADAPTER CLIP

1. Align the primary camera of the phone to the center of the circular hole of the adapter clip. Typically, the primary is a 1x wide-angle camera and the other camera(s) are secondary.

For the latest instructions, and for selecting primary camera for iPhone 13 Pro and later, visit www.carson.com/multicam




2. Attach the adapter clip to the top of the microscope and turn on camera. If viewing image is not centered or is not a complete round image, adjust adapter clip making sure it is centered over camera.

Note: Always hold the phone securely during use. If you release the phone while it is attached to the microscope it can fall and damage the phone.



3. Pinch screen to enlarge image. Take photos and videos to share with friends and family.

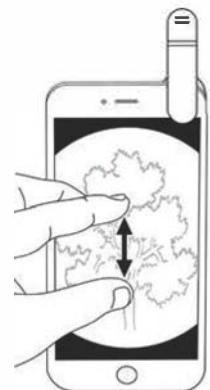

WARNING

1. To avoid permanent eye damage:
 - Do not look at the sun with the microscope.
 - Do not place the lighted microscope directly upon human eye.
2. Remove battery before storing microscope for extended periods.
3. The device may get warm after extended use. This is normal and should not be considered a defect.

Customer Service:

If you experience any difficulties, please contact us and we will be happy to help you.



✉ info@carson.com



✉ uksupport@carson.com



✉ eusupport@carson.com

For warranty information, visit www.carson.com/warranty

www.carson.com

Carson Optical 2070 5th Avenue, Ronkonkoma, NY 11779 | Tel: 631-963-5000